WDZX系列 >>>

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Intelligent microcomputer control system:
Using LED digital control system; After sterilization finished, sterilizer will automatic alarm and shut down.

Interlock device:
The control system can real time supervision the sterilizer chamber's temperature and pressure. Only under the safety temperature and pressure can open the sterilizer's cover. Make sure the operator's safety.

The structure of cover's opening:
Adopt turntable type rotary locking structure, make the lid open and lock safe, reliable, flexible and lightweight.

Preset program:
Can preset fixed program that sterilize solid body.

Memory storage system:
Store the last operation program so convenient next operation.

Feed water control:
Automatic water intake and water refill function.

Exhaust steam control :
After sterilization, exhaust steam valve should be opened (Digital display screen reminds exhaust steam) when sterilization time run out.

Interface check:
Provide temperature and pressure check interface; convenient to check multiple position temperature measurement.

Digital display screen:
Digital display screen and working indicator light display running state.

Safety devices:
Adopt the safety interlock structure of electrical and mechanical interaction to ensure that the interlock device locks the lid handle when there is pressure in the pot to avoid opening the lid when there is pressure in the pot.

Print function (add configuration) :
Print the current temperature and sterilization time.